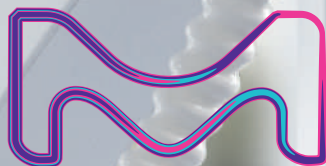
Milli-Q®

Lab Water

# Milli-Q® IQ 7000

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| 法規                                                   | 水質等級                                                        |
|------------------------------------------------------|-------------------------------------------------------------|
| European Pharmacopeia                                | Highly purified water                                       |
| U.S. Pharmacopeia                                    | Purified Water in bulk                                      |
| Japanese Pharmacopeia                                | Purified Water                                              |
| Chinese Pharmacopeia                                 | Purified Water                                              |
| ASTM® D1193                                          | Type I water                                                |
| ASTM® D5196                                          | Bio-Application Grade Water (with appropriate final filter) |
| ISO® 3696                                            | Grade 1 water                                               |
| Chinese National Standard                            | GB 6682 Grade 1 water                                       |
| JIS K 0557                                           | A4 water                                                    |
| Clinical and Laboratory Standards Institute® (CLSI®) | Clinical Laboratory Reagent Water (CLRW)                    |

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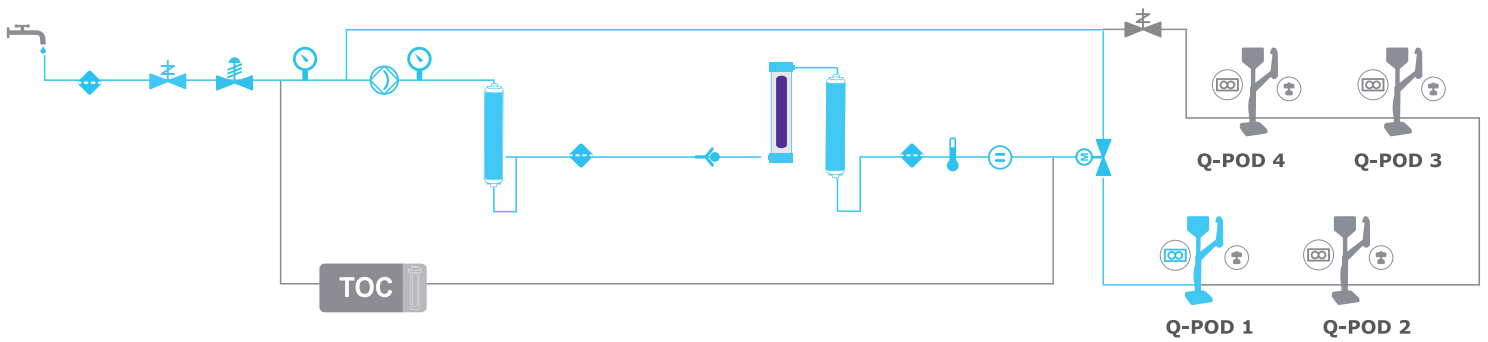
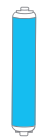
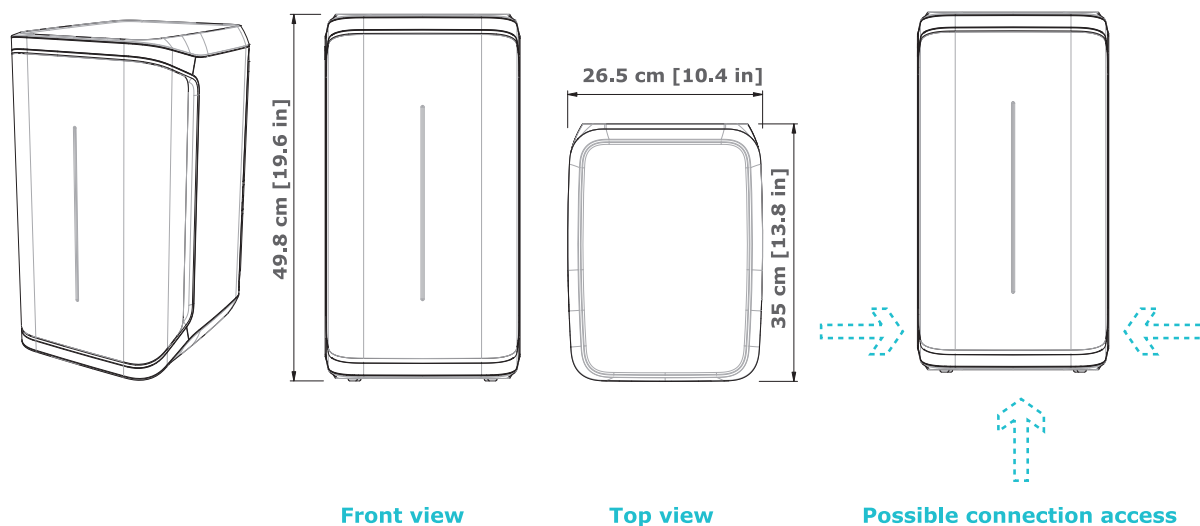
## Technical Appendix-2

### Tubing and port requirements

| Item                                                  | Description                                                      |
|-------------------------------------------------------|------------------------------------------------------------------|
| Feed water port                                       | PE tube dia 6x8 mm, maximum 5 m (16.4 ft) from feed              |
| Water to Q-POD® dispenser and back to production unit | PE tube dia 6x8 mm (inside connector sleeve), maximum length 5 m |
| Power entry                                           | Connection IEC 13                                                |
| ON/OFF switch                                         | Available on the unit                                            |
| Water sensor port                                     | Maximum 3.3 VDC                                                  |
| Tank level adapter port                               | Maximum 5 VDC                                                    |
| Ethernet port                                         | IEEE P802.3                                                      |

### Electrical connections and specifications

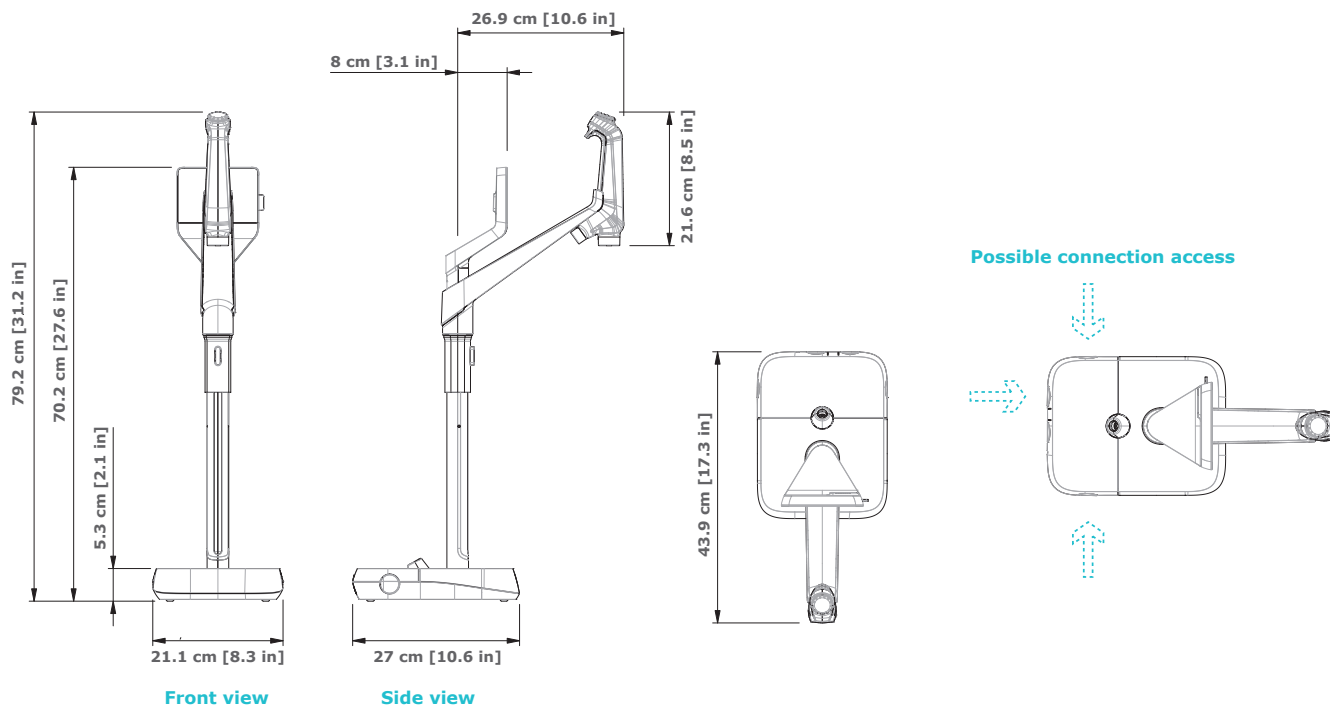
| Item                    | Description                              |
|-------------------------|------------------------------------------|
| Power source voltage    | 100 – 240 V ± 10%                        |
| Power frequency         | 50 – 60 Hz at ± 2Hz                      |
| Power used              | 350 VA                                   |
| Power cord              | Length 2.5 m (8.2 ft) Plug: IEC13 female |
| Operational temperature | 4 – 40 °C (39 – 104 °F)                  |
| Altitude                | 3000 m (9842 ft)                         |
| RFID frequency          | 13.56 MHz                                |

### Weight (Q-POD not included)

| Shipping weight | Dry weight          | Operating weight   |
|-----------------|---------------------|--------------------|
| 15 kg (33 lb)   | 12.46 kg (27.47 lb) | 16.06 kg (35.4 lb) |

# Q-POD® Dispenser Specifications

## Technical Appendix-3

### Tubeing and port requirements

| Item                                                                                           | Description                                                                                |
|------------------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------------|
| Dispenser tubing length                                                                        | 0.92 m (36.22 in)                                                                          |
| Distance from production unit to Q-POD® dispenser                                              | Maximum 5 m (16.4 ft)                                                                      |
| Distance between two Q-POD® dispensers<br>(Maximum four Q-POD® dispensers connected in series) | Maximum 5 m (16.4 ft)                                                                      |
| Q-POD® dispenser data connection with unit                                                     | Ethernet (cable inside connector sleeve)                                                   |
| Q-POD® electrical connection                                                                   | Q-POD® is powered by the production unit (24 VDC - 28 VDC) (cable inside connector sleeve) |
| RFID frequency                                                                                 | 13.56 MHz                                                                                  |
| Foot pedal port                                                                                | 3.3 V                                                                                      |

### Weight

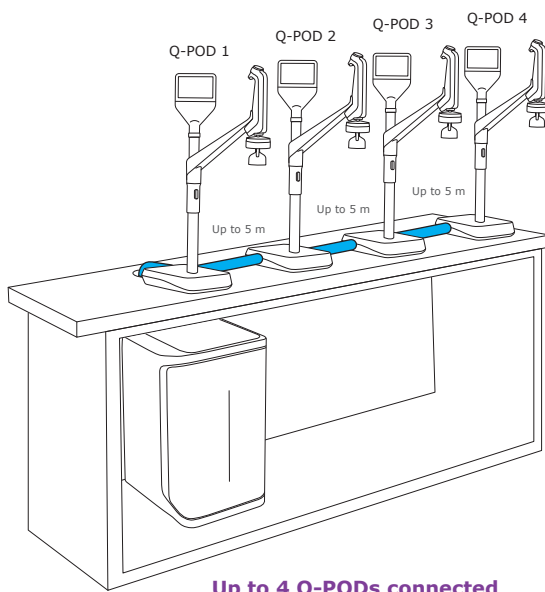
| Shipping weight  | Dry weight       | Operating weight  |
|------------------|------------------|-------------------|
| 7.9 kg (17.4 lb) | 5.5 kg (12.1 lb) | 5.64 kg (12.4 lb) |

### Screen description & functionalities


| Item                    | Description                                                                               |
|-------------------------|-------------------------------------------------------------------------------------------|
| Capacitive touch screen | Size: 5"<br>Resolution: 800 x 480                                                         |
| USB port                | USB 2.0 Highspeed standard                                                                |
| Speaker                 | Impedance: 8 Ω / Max output power: 0.5W                                                   |
| Display in 9 languages  | Chinese / English / French / German / Italian / Japanese / Portuguese / Russian / Spanish |

# Installation Options To Fit Your Requirements


## Technical Appendix-4


**Up to 4 Q-PODs connected**

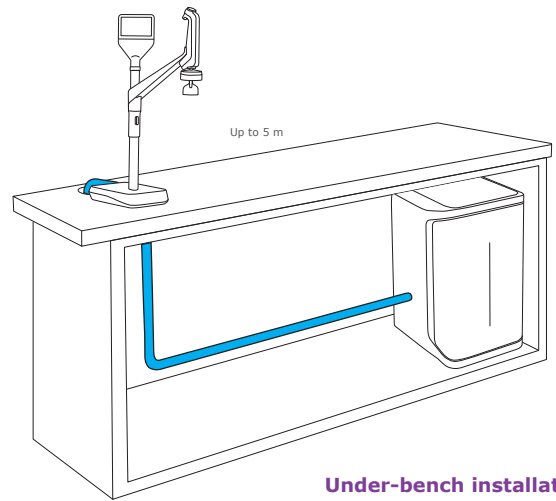**Under-bench installation**

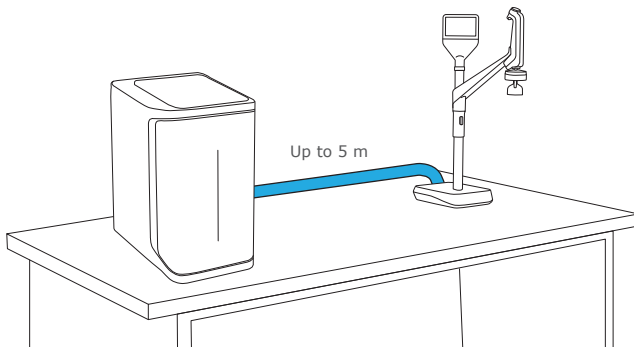**Bench-top installation**

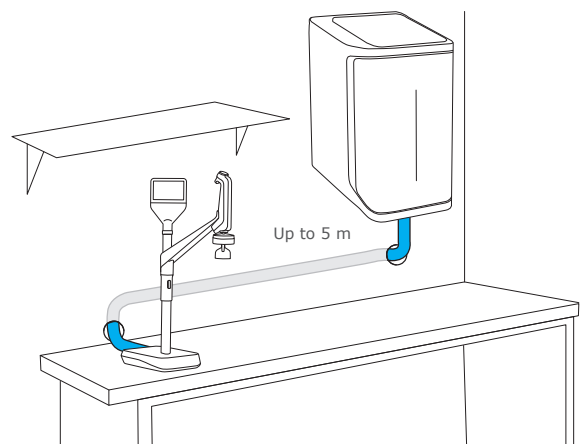**Wall-mounted installation**

# Water Specifications International Conformity

## Technical Appendix-5

### Feed water requirements

|                            |                                                                 |
|----------------------------|-----------------------------------------------------------------|
| Feed water quality         | Pretreated water using EDI, DI, RO or distillation technologies |
| Conductivity at 25 °C      | <100 µS/cm                                                      |
| Total Organic Carbon (TOC) | <50 ppb                                                         |
| Pressure range*            | Maximum 6 bar                                                   |
| Temperature                | 5 – 35 °C / 41 – 95 °F                                          |

\* For pressures between 0 and 0.1 bar, the system will operate, but product flow rate may be lower.

### Ultrapure, Type 1 water specifications<sup>(1)</sup> (from Q-POD dispenser)

|                                      |                                                               |
|--------------------------------------|---------------------------------------------------------------|
| Resistivity at 25 °C <sup>(2)</sup>  | 18.2 MΩ·cm                                                    |
| TOC                                  | ≤2 ppb (µg/L) <sup>(3)</sup> , typically ≤ 5 ppb (µg/L)       |
| Particles <sup>(4)</sup>             | No particles with size > 0.22 µm                              |
| Bacteria                             | <0.01 CFU/mL <sup>(4,5)</sup><br><0.005 CFU/mL <sup>(6)</sup> |
| Pyrogens (endotoxins) <sup>(5)</sup> | <0.001 EU/mL                                                  |
| RNases <sup>(5)</sup>                | <1 pg/mL                                                      |
| DNases <sup>(5)</sup>                | <5 pg/mL                                                      |
| Proteases <sup>(5)</sup>             | <0.15 µg/mL                                                   |
| Flow rate                            | 0.05 – 2 L/min                                                |

(1) These values are typical and may vary depending on the nature and concentration of contaminants in the feed water.

(2) Resistivity can also be displayed non-temperature-compensated as required by USP.

(3) In the appropriate operating conditions; otherwise typically ≤ 5 ppb.

(4) With Millipak® and Millipak® Gold.

(5) With Biopak®.

(6) With Millipak® Gold installed and used in a laminar flow hood.

## International regulation requirements

### Declaration of CE and cUL Conformity:

The Milli-Q® IQ 7000 system has been tested by an independent and accredited company for compliance with CE directives related to safety and

electromagnetic compatibility. The report can be consulted on demand. In addition, the Milli-Q IQ 7000 system is built using components and practices recommended by UL and has been cUL marked. The registration can be verified on the UL web site: [www.ul.com](http://www.ul.com)

**We also meet the regulatory requirements of the following organizations:**

# Ordering Information

## Technical Appendix-6

| Water purification system                  | Catalog number |
|--------------------------------------------|----------------|
| Milli-Q® IQ 7000 purification system       | ZIQ7000WW*     |
| Q-POD® remote dispenser with 2 m connector | ZIQP0D020      |
| Q-POD® remote dispenser with 5 m connector | ZIQP0D050      |

| Purification cartridges                           | Catalog number |
|---------------------------------------------------|----------------|
| IPAK Meta® & IPAK Quanta® polishing cartridge kit | IPAKKITA1      |
| Millipak® 0.22 µm filter                          | MPGP002A1      |
| Millipak® Gold 0.22 µm sterile filter             | MPGPG02A1      |
| Biopak® polisher                                  | CDUFBI0A1      |
| LC-Pak® polisher                                  | LCPAK00A1      |
| EDS-Pak® polisher                                 | EDSPAK0A1      |
| VOC-Pak® polisher                                 | V0CPAK0A1      |

| Accessories                                    | Catalog number |
|------------------------------------------------|----------------|
| Additional Q-POD® dispenser with 2 m connector | ZIQP0DX20      |
| Additional Q-POD® dispenser with 5 m connector | ZIQP0DX50      |
| Feed water conductivity cell                   | ZFC0NDCA1      |
| Foot pedal                                     | ZMQSFTSA1      |
| Wall mounting bracket                          | SYSTFIXA1      |
| Water sensor                                   | ZWATSENA1      |

| Services |
|----------|
|----------|

Our certified Field Service Engineers provide expert on-site professional support for the installation, validation, calibration and maintenance of your individual and total water solutions. To select the proper service plan for your water system, please contact your local service specialist to discuss and help you design the plan best adapted to your needs.

\* The end letters "WW" will be replaced by your country letters on your invoice to ensure you receive the appropriate accessory kit for your country, including the user manual in your local language and a power cord adapted to the local electrical network.

To order consumables online, please visit:  
[www.mymilliqconsumables.com](http://www.mymilliqconsumables.com)

For more information, please visit our website:  
[www.merckmillipore.com/milliq-iqsystem](http://www.merckmillipore.com/milliq-iqsystem)

For more information, please visit our website:  
[www.merckmillipore.com/labwater](http://www.merckmillipore.com/labwater)

Merck, MilliporeSigma, the vibrant M, Millipore, Milli-Q, Q-POD, Elix, Jetpore, ech<sub>2</sub>o, IPAK Meta, IPAK Quanta, IQnano, VOC-Pak, Millipak, Biopak, EDS-Pak, LC-Pak, and A10 are trademarks of Merck KGaA.

ISO is a registered trademark of the International Organization for Standardization.

Clinical and Laboratory Standards Institute and CLSI are registered trademarks of the Clinical and Laboratory Standards Institute, Inc.

ASTM is a registered trademark of the American Society for Testing and Materials.

All other trademarks are the property of their respective owners.

Lit. No. PB5475ENEU

© 2017 EMD Millipore Corporation, Billerica, MA, U.S.A. All rights reserved.